

## MestReNova (Mnova)

MestReNova (Mnova) is available to all University of Utah students and employees. Using Mnova requires downloading both the software and the main license file. Instructions for downloading and installing the software and license are provided below.

---

### Installation Instructions

#### 1. Download Mnova Software

- Our current license supports downloads up to revision 15.1.
- Versions are available for Mac, PC, and Linux.
- Download Mnova from this link: [Mnova Download](#).
- Select the version appropriate for your operating system.

#### 2. Download the License File

- Download the license file linked below:  
**Mnova License File**
- **PC Users:**
  - Right-click on the license file link and select "**Save Link As...**" from the context menu.
  - Choose a location to save the file. It is recommended to create a dedicated folder (e.g., Mnova\_License) for storing the .lic file. Alternatively, save the file to your **Downloads** folder.
- **Mac Users:**
  - Hold down the **Control** key and click on the license file link. From the menu that appears, select "**Download Linked File As...**" (or similar, depending on your browser).
  - Save the file in a dedicated folder (e.g., Mnova\_License) or in the default **Downloads** folder.
- **Important Note for All Users:**
  - Ensure the license file is saved with the .lic file extension intact. Do not rename or modify the file after downloading.
  - If your browser appends extra characters (e.g., .txt), manually rename the file to remove them, so it ends with .lic.
  - **A GlobalProtect VPN connection is required to validate the license so that the software can be used in most circumstances. ##add link to VPN instructions page here**

#### 3. Install Mnova

- Begin by installing the Mnova software.
- Once the installation is complete, start Mnova and follow the on-screen instructions to install the license file.

#### 4. Optional Plugins

- Optional plugins can be installed using the Mnova Registration Wizard. Links to the plugin license files are provided below:
  - **[Mnova qNMR License File]**
  - **[Mnova Reaction Monitor License File]**